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cDNA synthesis kit (CN: KPG-cDNA50)

A short explanation about the cDNA synthesis kit

Considering that RNA is unstable at high temperatures, it must be converted to DNA for use in some techniques such as PCR. The cDNA synthesis kit of Karmania Pars Gene Company contains all the necessary components to convert total RNA or mRNA into cDNA. The current kit has been designed and produced using Reverse Transcriptase enzyme of Karmania Pars Gene Company. This enzyme is designed and expressed in such a way that it is resistant to temperature and lacks RNase H properties. On the other hand, it has an optimal performance compared to the enzymes available in the Iranian market. The next feature of this kit is ease of use for the consumer and it can be done with minimal sampling.

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Procedure

For cDNA synthesis, add the following to the RNase/DNase free microtube.

Template RNA	5 μL recommended
Oligo dt or Random hexamer	1 μL
cDNA Master mix	14 μL
Total	Up to 20 μL

RNA between 5 ng and 5 μ g can be used. After adding the above, first mix a little and then run the following temperature program:

Temperature	Time
25°C	15 minutes
37°C	15 minutes
45°C	30 minutes
55°C	15 minutes
80°C	10 minutes

To check the accuracy of the kit during PCR, use the beta-actin primer in the kit as a positive control. To perform PCR, run 1 microliter of primer in the PCR mixture with the following temperature program:

Cycle	Temperature	Time
1	95°C	5 minutes
35	95°C	20 Seconds
	60°C	25 Seconds

The temperature program is based on the use of Master Biosystem PCR master mix. If you use Mastermix Cybergreen from other companies, use 58°C for the annealing temperature and follow the temperature program of Mastermix. Our recommended program:

Cycle	Temperature	Time
1	95°C	2 minutes
35	95°C	20 Seconds
	58°C	25 Seconds
	72°C	30 Seconds

Safety:

The solutions used in the kit are not harmful for humans, however, avoid direct contact with solutions.

Contact us

Block 269, Food and Drug area, Rafsanjan Special Economic Zone, Rafsanjan-Iran

Tel: +983434208024-25, +989132926113, +989135025983

E-Mail: Karmaniaparsgene@gmail.com

Website: www.karmaniaparsgene.ir

Samples:

You must use purified RNA for cDNA synthesis. Our suggestion is using KPG RNA Extraction kits

Other PCR-related products of Karmania Pars Gene Company:

Name	Catalog Number
PCR Master Mix (2 vials)	(CN: KPG-MM)
Agarose	(CN:KPG-AG)
50 bp Ladder	(CN:KPG-La50)
100 bp Ladder	(CN:KPG- La100)
1kb Ladder	(CN:KPG- La1k)
Thermoresistance RT	(CN:KPG-trt)

The list of products in this area is being updated, Please check our website.

Other products of KPG Company:

ELISA kits to measure cytokines
Oxidant and Antioxidant Kits
Real-Time PCR Microtubes
RNA, DNA, microRNA and Plasmid Extraction kits
Real-Time PCR Pathogens Testing Kits
DNA, RNA Spin Column
Collection Tube
Filter head Sampler Tips
Stem, Cancer and primary Cell Lines

Visit our website for the latest updated list of products.

Website: www.karmaniaparsgene.ir